



PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit

Catalog No.	Quantity	Total Purifications
12955-4	4 Preps	384
12955-12	12 Preps	1152

Instruction Manual

Inhibitor Removal Technology[®] (IRT) is a registered trademark of MO BIO Laboratories, Inc. and is covered by US patent protection as well as international patents pending.



Please recycle

Version: 01062012

Technical Information: Toll free 1-800-606-6246, or 1-760-929-9911 Email: technical@mobio.com Website: www.mobio.com



Table of Contents

Introduction	3
Protocol Overview	3
Equipment Required	4
Kit Contents & Storage	4
Precautions & Warnings	4
Protocols:	
Experienced User Combination Vacuum & Centrifugation Protocol	5
Detailed Combination Protocol (Describes what is happening at each step)	7
Experienced User Centrifugation Protocol	11
Detailed Centrifugation Protocol (Describes what is happening at each step)	13
Hints & Troubleshooting Guide	17
Additional Applications and References	19
Contact Information	21
Other Quality Products Available	22



Introduction

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit can be used for isolating genomic DNA from up to 0.25 grams of soil.

Two protocols are included for use with this product. The first protocol involves a combination of vacuum and centrifugation steps. The second protocol uses centrifugation exclusively. This kit requires the use of a specialized plate shaker for the process of shaking beads in deep well plates which aid in cell lysis. We recommend the Retsch plate shaker (MO BIO Catalog# 11996).

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit provides researchers with a novel and proprietary high throughput method for isolating genomic DNA from environmental samples. The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit distinguishes itself from MO BIO Laboratories' UltraClean[®]-htp 96 Well Soil DNA Isolation Kit with a patented Inhibitor Removal Technology[®] (IRT) procedure that is much more efficient at removing humic substances that inhibit PCR. This new procedure is effective at removing PCR inhibitors from even the most difficult soil types allowing for more successful PCR amplification of DNA.

Tested sample types include compost, sediment, and manure among many others. Isolated DNA was successfully used as template DNA to amplify members of a wide range of microbial groups in soils. These include bacteria (gram positive, gram negative, and spore formers), actinomycetes, and fungi.

Protocol Overview

Environmental samples are added to a 96 well bead beating plate for rapid and thorough homogenization. Cell lysis occurs by a combination of mechanical and chemical methods. Humics are removed by a specialized precipitation process. Total genomic DNA is captured on a 96 well silica membrane in a spin column plate format. DNA is then washed and eluted from the membrane. The eluted DNA is ready for PCR analysis and other downstream applications.

NOTE: The estimated time from start to finish required to process two 96 well plates for this protocol is approximately 8 hours. Stopping points at appropriate steps are mentioned in the protocol. The majority of the time is for weighing and loading the soil samples into the 96 well plates.

This kit is for research purposes only. Not for diagnostic use.

Other Related Products	Catalog No.	Quantity
96 Well Plate Shaker	11996	1 unit (120 V)
	11996-220	1 unit (220 V)
Vacuum Pump	11998	1 unit (120 V)
	11998-220	1 unit (220 V)
Plate Adapter Set	11999	1 set



Equipment Required

Centrifuge capable of handling two 96 Well blocks (13 cm x 8.5 cm x 6.0 cm) at 4500 x g

Note: If you have a centrifuge with a maximum speed less than 4500 x g see the Hints and Troubleshooting Guide.

Multi-channel Pipettor (volumes of 50 µl - 650 µl)

Mechanical Shaker for 96 Well Blocks and Plate Adapters (MO BIO Catalog# 11996 and 11999)

Vortex-Genie[®] 2 Vortex with 3 inch platform (MO BIO Catalog# 13111-V or 13111-V-220)

Reagents Required but not Included

100% Ethanol

Optional Material/Equipment

Reagent reservoirs

Vacuum Pump (MO BIO Catalog# 11998)

Vacuum Manifold (MO BIO recommends Axygens' AxyVac Vacuum Manifold Catalog# AP-VM)

Kit Contents

Component	Kit Catalog # 12955-4		Kit Catalog # 12955-12	
	Catalog #	Amount	Catalog #	Amount
PowerSoil [®] -htp Bead Plate (w/Square Well Mat)	12955-4-BP	4	12955-12-BP	12
PowerSoil [®] -htp Bead Solution	12955-4-BS	340 ml	12955-12-BS	1020 ml
PowerSoil [®] -htp Solution C1	12955-4-1	45 ml	12955-12-1	135 ml
PowerSoil [®] -htp Solution C2	12955-4-2	128 ml	12955-12-2	385 ml
PowerSoil [®] -htp Solution C3	12955-4-3	106 ml	12955-12-3	320 ml
PowerSoil [®] -htp Solution C4	12955-4-4	550 ml	12955-12-4	2 x 825 ml
PowerSoil [®] -htp Solution C5-D Before first use, Solution C5-D must be prepared. Please see step 1 of protocol.	12955-4-5D	120 ml	12955-12-5D	360 ml
PowerSoil [®] -htp Solution C6	12955-4-6	66 ml	12955-12-6	200 ml
PowerSoil [®] -htp Spin Plates	12955-4-SP	4	12955-12-SP	12
PowerSoil [®] -htp 2 ml Collection Plates	12955-4-2CP	8	12955-12-2CP	24
PowerSoil [®] -htp 1 ml Collection Plates	12955-4-1CP	16	12955-12-1CP	48
PowerSoil [®] -htp 0.5 ml Collection Plates	12955-4-0.5CP	4	12955-12-0.5CP	12
PowerSoil [®] -htp Microplates	12955-4-MP	4	12955-12-MP	12
PowerSoil [®] -htp Centrifuge Tape	12955-4-CT	24	12955-12-CT	72
PowerSoil [®] -htp Sealing Tape	12955-4-ST	16	12955-12-ST	48
PowerSoil [®] -htp Elution Sealing Mats	12955-4-ESM	4	12955-12-ESM	12

Kit Storage

Kit reagents and components should be stored at room temperature (15-30°C).

Precautions

Please wear gloves when using this product. Avoid all skin contact with kit reagents. In case of contact, wash thoroughly with water. Do not ingest. See Material Safety Data Sheets for emergency procedures in case of accidental ingestion or contact. All MSDS information is available upon request (760-929-9911) or at www.mobio.com. Reagents labeled flammable should be kept away from open flames and sparks.

WARNING: Solution C5-D is flammable after addition of ethanol.

Technical Information: Toll free 1-800-606-6246, or 1-760-929-9911 Email: technical@mobio.com Website: www.mobio.com

Experienced User Combination Vacuum and Centrifugation Protocol

Please wear gloves at all times

- BEFORE THE FIRST USE ONLY, Solution C5-D must be prepared. Add an equal volume of 100% Ethanol to Solution C5-D (for the 4 prep kit = 120 ml, or for the 12 prep kit = 360 ml). Mix well. Put a check mark in the “ethanol added” box on the bottle cap label.**
- Remove the Square Well Mat from the PowerSoil[®]-htp Bead Plate and set aside. Add 0.25 grams of soil sample.
Note: This is an appropriate stopping point and you can store the PowerSoil[®]-htp Bead Plate at 4°C covered with the Square Well Mat.
This is the most time consuming step of the protocol. Care must be taken to avoid cross contamination between sample wells.
- Add 750 µl of PowerSoil[®]-htp Bead Solution to the wells of the PowerSoil[®]-htp Bead Plate.
- Check Solution C1.** If Solution C1 has precipitated, heat the solution at 60°C until the precipitate has dissolved. Mix gently before using.
Note: *Solution C1 contains SDS. If it gets cold, it will precipitate. Heating at 60°C will dissolve the SDS. Solution C1 can be used while it is still warm.*
- Add 60 µl of Solution C1. Secure the Square Well Mat (from step 2) tightly to the PowerSoil[®]-htp Bead Plate.
Note: *A proper seal of the mat is critical to prevent loss of sample and leakage that might cause damage to your shaker.*
- Place PowerSoil[®]-htp Bead Plate with Square Well Mat securely fastened in the 96 Well Plate Shaker (MO BIO Catalog# 11996).
Note: The final order of assembly is; Aluminum Plate, Square Well Mat, 96 Well PowerSoil[®]-htp Bead Plate, and second Aluminum Plate.
- Shake at speed 20 for 10 minutes, remove plates and re-orient them so that the side closest to the machine body is now furthest from the machine body and shake again at speed 20 for 10 minutes.
- Centrifuge at room temperature for 6 minutes at 4500 x g.
- Remove the Bead Plate from the centrifuge. Carefully and without splashing remove and discard the Square Well Mat and transfer the supernatant to a clean 1 ml Collection Plate.
Note: The supernatant may still contain some soil particles.
- Add 250 µl of Solution C2 to each well and apply Sealing Tape to 1 ml Collection Plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge the 1 ml Collection Plate at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.
- Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.
- Apply new Sealing Tape to the 1 ml Collection Plate and centrifuge at room temperature for 6 minutes at 4500 x g. Transfer entire volume of supernatant to a new 1 ml Collection Plate.
- Add 200 µl of Solution C3 and apply Sealing Tape to the 1 ml Collection Plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.
- Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.
- Apply Sealing Tape to 1 ml Collection Plate. Centrifuge the 1 ml Collection Plate at room temperature for 6 minutes at 4500 x g.
- Transfer no more than **650 µl** of supernatant to a 2 ml Collection Plate avoiding any residual pellet.
- Add 650 µl of Solution C4 to each well of the 2 ml Collection Plate.
- Add a second 650 µl of Solution C4 to each well of the 2 ml Collection Plate.
Note: It is safe to stop the protocol at this step and store the samples covered with Sealing Tape, at 4°C.



19. Remove the top portion of the vacuum manifold and place a new 2 ml Collection Plate in the bottom of the vacuum manifold.
Note: If you are using the Axygen AP-VM, use the Waste Reservoir provided with the manifold in place of the 2 ml Collection Plate.
20. Replace the top of the manifold and place a Spin Plate on top of the manifold and turn the vacuum pump on.
21. Test that you have a good seal with the manifold and the Spin Plate by gently lifting up on the Spin Plate. You should be able to lift the entire unit without the Spin Plate separating from the manifold.
22. Pipet samples from step 18 “up and down” to mix.
23. Load approximately 650 μ l into the wells of each row of the Spin Plate until all the samples have been processed. Allow the liquid to pass through the wells.
24. Once the liquid has passed through the wells, continue loading 650 μ l volumes into the wells of each row of the Spin Plate until the entire volume of each sample is processed. Turn off the vacuum.
25. Remove the Spin Plate from the manifold and set aside, discard the flow through from the 2 ml Collection Plate or Waste Reservoir in the bottom of the manifold and then place it back in the manifold.
26. Reassemble the manifold with the Spin Plate back on top of the manifold and turn the vacuum on. Test the seal by lifting gently.
27. **Confirm that ethanol has been added to Solution C5-D (see step 1).** Add 500 μ l of Solution C5-D to each well of the Spin Plate. After the entire volume of Solution C5-D has passed through the wells of the Spin Plate, turn off the vacuum.
28. Apply Centrifuge Tape to the Spin Plate to cover the wells.
29. Place a 0.5 ml Collection Plate under the Spin Plate. Centrifuge at room temperature for 6 minutes at 4500 x *g*.
30. Carefully place Spin Plate onto a Microplate. Remove Centrifuge Tape and discard.
31. Allow to air dry for 10 minutes at room temperature.
32. Add 100 μ l of Solution C6 to the center of each well of the Spin Plate and apply Centrifuge Tape.
33. Centrifuge at room temperature for 3 minutes at 4500 x *g*. Remove Centrifuge Tape and discard.
34. Cover wells of Microplate with the Elution Sealing Mat provided. DNA is now ready for any downstream application. No further steps are required.

Prolonged storage at 4°C will result in the evaporation of eluted DNA. We recommend storing DNA frozen (-20°C or -80°C). Solution C6 does not contain EDTA. To concentrate the DNA see the Hints and Troubleshooting Guide.

Thank you for choosing the PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit.



Detailed Combination Vacuum and Centrifugation Protocol (Describes what is happening at each step)

Please wear gloves at all times

1. **BEFORE THE FIRST USE ONLY, Solution C5-D must be prepared. Add an equal volume of 100% Ethanol to Solution C5-D (for the 4 prep kit = 120 ml, or for the 12 prep kit = 360 ml). Mix well. Put a check mark in the "ethanol added" box on the bottle cap label.**

2. Remove the Square Well Mat from the PowerSoil[®]-htp Bead Plate and set aside. Add 0.25 grams of soil sample.

Note: This is an appropriate stopping point and you can store the PowerSoil[®]-htp Bead Plate at 4°C covered with the Square Well Mat.

What's happening: This is the most time consuming step of the protocol. Care must be taken to avoid cross contamination between sample wells.

3. Add 750 µl of PowerSoil[®]-htp Bead Solution to the wells of the PowerSoil[®]-htp Bead Plate.

What's happening: The soil sample has now been loaded into the PowerSoil[®]-htp Bead Plate. This is the first part of the lysis procedure. The PowerSoil[®]-htp Bead Solution is a buffer that will disperse the soil particles and begin to dissolve humic acids.

4. **Check Solution C1.** If Solution C1 has precipitated, heat the solution at 60°C until the precipitate has dissolved.

What's happening: Solution C1 contains SDS. If it gets cold, it will precipitate. Heating at 60°C will dissolve the SDS. Solution C1 can be used while it is still warm.

5. Add 60 µl of Solution C1. Secure the Square Well Mat (from step 2) tightly to the PowerSoil[®]-htp Bead Plate.

What's happening: Solution C1 contains SDS. This is a detergent that aids in cell lysis. The detergent breaks down fatty acids and lipids associated with the cell membrane of several organisms.

6. Place PowerSoil[®]-htp Bead Plate with Square Well Mat securely fastened (from step 2), to the 96 Well Plate Shaker (MO BIO Catalog# 11996).

Note: The final order of assembly is; Aluminum Plate, Square Well Mat, 96 Well PowerSoil[®]-htp Bead Plate, and second Aluminum Plate.

What's happening: Mechanical lysis is introduced at this step. The protocol uses a combination of mechanical and chemical lysis. By randomly shaking the beads, they collide with one another and with microbial cells causing them to break open.

7. Shake at speed 20 for 10 minutes, remove plates and re-orient them so that the side closest to the machine body is now furthest from the machine body and shake again at speed 20 for 10 minutes.

8. Centrifuge at room temperature for 6 minutes at 4500 x g.

What's happening: Particulates including cell debris, soil, beads, and humic acids, will form a pellet at this point. DNA is in the liquid supernatant.

9. Remove the Bead Plate from the centrifuge. Carefully and without splashing, remove and discard the Square Well Mat and transfer the supernatant to a clean 1 ml Collection Plate.

Note: The supernatant may still contain some soil particles.

10. Add 250 μ l of Solution C2 to each well and apply Sealing Tape to 1 ml Collection Plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge the 1 ml Collection Plate at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.

What's happening: Solution C2 contains a reagent to precipitate non-DNA organic and inorganic material including humic substances, cell debris, and proteins. It is important to remove contaminating organic and inorganic matter that may reduce DNA purity and inhibit downstream DNA applications.

11. Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate. Apply Sealing Tape to the 1 ml Collection Plate. Centrifuge the 1 ml Collection Plate at room temperature for 6 minutes at 4500 x g. Transfer entire volume of supernatant to another new 1 ml Collection Plate.

12. Add 200 μ l of Solution C3 and apply Sealing Tape to the 1 ml Collection Plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.

What's happening: Solution C3 is a second reagent to precipitate additional non-DNA organic and inorganic material including humic substances, cell debris, and proteins. It is important to remove contaminating organic and inorganic matter that may reduce DNA purity and inhibit downstream DNA applications.

13. Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.

What's happening: The pellet at this point contains residues of humic substances, cell debris, and proteins. For the best DNA yield and quality, avoid transferring any of the pellet.

14. Apply Sealing Tape to the 1 ml Collection Plate. Centrifuge the 1 ml Collection Plate at room temperature for 6 minutes at 4500 x g.

15. Transfer no more than **650 μ l** of supernatant to a 2 ml Collection Plate avoiding any residual pellet.

16. Add 650 μ l of Solution C4 to each well of the 2 ml Collection Plate.

What's happening: Solution C4 is a high concentration salt solution. Since DNA binds tightly to silica at high salt concentrations, this solution will adjust the salt concentration to allow binding of DNA, but not non-DNA organic and inorganic material that may still be present at low levels, to the Spin Plate.

17. Add a second 650 μ l of Solution C4 to each well of the 2 ml Collection Plate.

Note: It is safe to stop the protocol at this step and store the samples covered with Sealing Tape, at 4°C.

18. Remove the top portion of the vacuum manifold and place a new 2 ml Collection Plate in the bottom of the vacuum manifold.

Note: If you are using the Axygen AP-VM, use the Waste Reservoir provided with the manifold in place of the 2 ml Collection Plate.



19. Replace the top of the manifold and place a Spin Plate on top of the manifold and turn the vacuum pump on.
20. Test that you have a good seal with the manifold and the Spin Plate by gently lifting up on the Spin Plate. You should be able to lift the entire unit without the Spin Plate separating from the manifold.
21. Pipet samples from step 17 “up and down” to mix.
22. Load approximately 650 μ l into the wells of each row of the Spin Plate until all the samples have been processed. Allow the liquid to pass through the wells.
23. Once the liquid has passed through the wells continue loading 650 μ l volumes into the wells of each row of the Spin Plate until the entire volume of each sample is processed. Turn off the vacuum.

What's happening: DNA is selectively bound to the silica membrane in the Spin Plate device in the high salt solution. Contaminants pass through the filter membrane, leaving only DNA bound to the membrane.

24. Remove the Spin Plate from the manifold and set aside, discard the flow through from the 2 ml Collection Plate or Waste Reservoir in the bottom of the manifold and then place it back in the manifold.
25. Reassemble the manifold with the Spin Plate back on top of the manifold and turn the vacuum on. Test the seal by lifting gently.
26. **Confirm that ethanol has been added to Solution C5-D (see step 1).** Add 500 μ l of Solution C5-D to each well of the Spin Plate. After the entire volume of Solution C5-D has passed through the wells of the Spin Plate, turn off the vacuum.

What's happening: Solution C5-D is an ethanol based wash solution used to further clean the DNA that is bound to the silica membrane in the Spin Plate. This wash solution removes residues of salt, humic acid, and other contaminants while allowing the DNA to stay bound to the silica membrane. This flow through is just waste containing ethanol wash solution and contaminants that did not bind to the silica Spin Plate membrane.

Note: You can wash more than one time to further clean DNA if desired. In some cases where soils have very high humic acid content, it will be beneficial to repeat this wash step. There is 10% extra Solution C5-D in the bottle for this purpose. Solution C5-D is also sold separately (MO BIO Catalog# 12955-4-5D & 12955-12-5D).

27. Apply Centrifuge Tape to the Spin Plate to cover the wells.
28. Place a 0.5 ml Collection Plate under the Spin Plate. Centrifuge at room temperature for 6 minutes at 4500 x g.

What's happening: This step removes residual Solution C5-D (ethanol wash solution). It is critical to remove all traces of wash solution because it can interfere with many downstream DNA applications.

29. Carefully place Spin Plate onto a Microplate. Remove Centrifuge Tape and discard. Allow to air dry for 10 minutes at room temperature.

What's happening: This step removes residual Solution C5-D (ethanol wash solution). It is critical to remove all traces of wash solution because it can interfere with many downstream DNA applications.

Technical Information: Toll free 1-800-606-6246, or 1-760-929-9911 Email: technical@mobio.com Website: www.mobio.com



30. Add 100 µl of Solution C6 to the center of each well of the Spin Plate and apply Centrifuge Tape.

What's happening: Placing the Solution C6 (sterile elution buffer) in the center of the small white membrane will make sure the entire membrane is wetted. This will result in more efficient and complete release of the DNA from the Spin Plate membrane.

31. Centrifuge at room temperature for 3 minutes at 4500 x g. Remove Centrifuge Tape and discard.

What's happening: As the Solution C6 (elution buffer) passes through the silica membrane, DNA that was bound in the presence of high salt is selectively released by Solution C6 (10 mM Tris pH 8.0) which is a low salt solution.

32. Cover wells of Microplate with the Elution Sealing Mat provided. DNA is now ready for any downstream application. No further steps are required.

Prolonged storage at 4°C will result in the evaporation of eluted DNA. We recommend storing DNA frozen (-20°C or -80°C). Solution C6 does not contain EDTA. To concentrate the DNA see the Hints and Troubleshooting Guide.

Thank you for choosing the PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit.



Experienced User Centrifugation Protocol

Please wear gloves at all times

- BEFORE THE FIRST USE ONLY, Solution C5-D must be prepared. Add an equal amount of 100% Ethanol to Solution C5-D (for the 4 prep kit = 120 ml, or for the 12 prep kit = 360 ml). Mix well. Put a check mark in the “ethanol added” box on the bottle cap label.**
- Remove the Square Well Mat from the PowerSoil[®]-htp Bead Plate and set aside. Add 0.25 grams of soil sample.
Note: This is an appropriate stopping point and you can store the PowerSoil[®]-htp Bead Plate at 4°C covered with the Square Well Mat.
This is the most time consuming step of the protocol. Care must be taken to avoid cross contamination between sample wells.
- Add 750 µl of PowerSoil[®]-htp Bead Solution to the wells of the PowerSoil[®]-htp Bead Plate.
- Check Solution C1.** If Solution C1 has precipitated, heat solution at 60°C until the precipitate has dissolved.
Note: *Solution C1 contains SDS. If it gets cold, it will precipitate. Heating at 60°C will dissolve the SDS. Solution C1 can be used while it is still warm.*
- Add 60 µl of Solution C1. Secure the Square Well Mat (from step 2) tightly to the plate.
- Place PowerSoil[®]-htp Bead Plate with Mat securely fastened to the 96 Well Plate Shaker (MO BIO Catalog# 11996).
Note: The final order of assembly is; Aluminum Plate, Square Well Mat, 96 Well PowerSoil[®]-htp Plate, and second Aluminum Plate.
- Shake at speed 20 for 10 minutes, remove plates and re-orient them so that the side closest to the machine body is now furthest from the machine body and shake again at speed 20 for 10 minutes.
- Centrifuge at room temperature for 6 minutes at 4500 x g.
- Remove and discard the Square Well Mat. Transfer the supernatant to a clean 1 ml Collection Plate.
Note: The supernatant may still contain some soil particles.
- Add 250 µl of Solution C2 and apply Sealing Tape to plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge the plate at room temperature for 17 minutes at 4500 x g. Remove and discard Sealing Tape.
- Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.
- Apply Sealing Tape to plate. Centrifuge the plate again at room temperature for 6 minutes at 4500 x g. Transfer entire volume of supernatant to another new 1 ml Collection Plate.
- Add 200 µl of Solution C3 and apply Sealing Tape to plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.
- Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.
- Apply Sealing Tape to plate. Centrifuge the plate again at room temperature for 6 minutes at 4500 x g.
- Transfer no more than **650 µl** of supernatant to a 2 ml Collection Plate avoiding any residual pellet.
- Add 650 µl of Solution C4 to each well of the plate.
- Add a second 650 µl of Solution C4 to each well of the plate.
Note: It is safe to stop the protocol at this step and store the samples covered with Sealing Tape at 4°C.
- Pipet samples “up and down” to mix.
- Place Spin Plate onto a new 0.5 ml Collection Plate.
- Load approximately 650 µl into each well of the Spin Plate and apply Centrifuge Tape.
- Centrifuge at room temperature for 3 minutes at 4500 x g. Discard the flow through and place the Spin Plate back on the same 0.5 ml Collection Plate. Discard the Centrifuge Tape.
- Repeat steps 20-21 until all the supernatant has been processed. Discard the final flow through.
- Place the Spin Plate back on the same 0.5 ml Collection Plate.

Technical Information: Toll free 1-800-606-6246, or 1-760-929-9911 Email: technical@mobio.com Website: www.mobio.com



25. **Confirm that ethanol has been added to Solution C5-D (see step 1).** Add 500 µl of Solution C5-D to each well of the Spin Plate. Apply Centrifuge Tape to the Spin Plate.
26. Centrifuge at room temperature for 3 minutes at 4500 x *g*. Discard the flow through and place the same 0.5 ml Collection Plate beneath the Spin Plate.
27. Centrifuge again at room temperature for 5 minutes at 4500 x *g*. Discard the flow through.
28. Carefully place the Spin Plate onto a Microplate. Remove Centrifuge Tape and discard.
29. Allow to air dry for 10 minutes at room temperature.
30. Add 100 µl of Solution C6 to the center of each well of the Spin Plate. Apply Centrifuge Tape.
31. Centrifuge at room temperature for 3 minutes at 4500 x *g*. Remove Centrifuge Tape and discard.
32. Cover wells of Microplate with the Elution Sealing Mat provided. DNA is now ready for any downstream application. No further steps are required.

Prolonged storage at 4°C will result in the evaporation of eluted DNA. We recommend storing DNA frozen (-20°C or -80°C). Solution C6 does not contain EDTA. To concentrate the DNA see the Hints and Troubleshooting Guide.

Thank you for choosing the PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit.

Detailed User Centrifugation Protocol (Describes what is happening at each step)

Please wear gloves at all times

1. **BEFORE THE FIRST USE ONLY, Solution C5-D must be prepared. Add an equal amount of 100% Ethanol to Solution C5-D (for the 4 prep kit = 120 ml, or for the 12 prep kit = 360 ml). Mix well. Put a check mark in the “ethanol added” box on the bottle cap label.**

2. Remove the Square Well Mat from the PowerSoil[®]-htp Bead Plate and set aside. Add 0.25 grams of soil sample.

Note: This is an appropriate stopping point and you can store the PowerSoil[®]-htp Bead Plate at 4°C covered with the Square Well Mat.

What's happening: This is the most time consuming step of the protocol. Care must be taken to avoid cross contamination between sample wells.

3. Add 750 µl of PowerSoil[®]-htp Bead Solution to the wells of the PowerSoil[®]-htp Bead Plate.

What's happening: The soil sample has now been loaded into the PowerSoil[®]-htp Bead Plate. This is the first part of the lysis procedure. The PowerSoil[®]-htp Bead Solution is a buffer that will disperse the soil particles.

4. **Check Solution C1.** If Solution C1 has precipitated, heat solution at 60°C until precipitate has dissolved.

What's happening: Solution C1 contains SDS. If it gets cold, it will precipitate. Heating at 60°C will dissolve the SDS. Solution C1 can be used while it is still warm.

5. Add 60 µl of Solution C1. Secure the Square Well Mat (from step 2) tightly to the plate.

What's happening: Solution C1 contains SDS. This is a detergent that aids in cell lysis. The detergent breaks down fatty acids and lipids associated with the cell membrane of microorganisms.

6. Place PowerSoil[®]-htp Bead Plate with Mat securely fastened, to the 96 Well Plate Shaker (MO BIO Catalog# 11996).

Note: The final order of assembly is; Aluminum Plate, Square Well Mat, 96 Well PowerSoil[®]-htp Plate, and second Aluminum Plate.

What's happening: Mechanical lysis is introduced at this step. The protocol uses a combination of mechanical and chemical lysis. By randomly shaking the beads, they collide with one another and with microbial cells causing them to break open.

7. Shake at speed 20 for 10 minutes, remove plates and re-orient them so that the side closest to the machine body is now furthest from the machine body and shake again at speed 20 for 10 minutes.

8. Centrifuge at room temperature for 6 minutes at 4500 x g.

What's happening: Particulates including cell debris, soil, beads, and humic acids, will form a pellet at this point. DNA is in the liquid supernatant.

9. Remove and discard the Square Well Mat. Transfer the supernatant to a clean 1 ml Collection Plate.

Note: The supernatant may still contain some soil particles.

10. Add 250 μ l of Solution C2 and apply Sealing Tape to plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge the plate at room temperature for 17 minutes at 4500 x g. Remove and discard Sealing Tape.

What's happening: Solution C2 contains a reagent to precipitate non-DNA organic and inorganic material including humic substances, cell debris, and proteins. It is important to remove contaminating organic and inorganic matter that may reduce DNA purity and inhibit downstream DNA applications.

11. Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.
12. Apply Sealing Tape to plate. Centrifuge the plate again at room temperature for 6 minutes at 4500 x g. Transfer entire volume of supernatant to another new 1 ml Collection Plate.
13. Add 200 μ l of Solution C3 and apply Sealing Tape to plate. Vortex for 5 seconds and incubate at 4°C for 10 minutes. Centrifuge at room temperature for 6 minutes at 4500 x g. Remove and discard Sealing Tape.

What's happening: Solution C3 is a second reagent to precipitate additional non-DNA organic and inorganic material including humic substances, cell debris, and proteins. It is important to remove contaminating organic and inorganic matter that may reduce DNA purity and inhibit downstream DNA applications.

14. Avoiding the pellet, transfer entire volume of supernatant to a new 1 ml Collection Plate.

What's happening: The pellet at this point contains residues of humic substances, cell debris, and proteins. For best DNA yield and quality, avoid transferring any of the pellet.

15. Apply Sealing Tape to plate. Centrifuge the plate again at room temperature for 6 minutes at 4500 x g.
16. Transfer no more than **650 μ l** of supernatant to a 2 ml Collection Plate avoiding any residual pellet.
17. Add 650 μ l of Solution C4 to each well of the plate.

What's happening: Solution C4 is a high concentration salt solution. Since DNA binds tightly to silica at high salt concentrations, this solution will adjust the salt concentrations to allow binding of DNA, but not non-DNA organic and inorganic material that may still be present at low levels, to the Spin Plate.

18. Add a second 650 μ l of Solution C4 to each well of the plate.

Note: It is safe to stop the protocol at this step and store the samples covered with Sealing Tape, at 4°C.

19. Pipet samples "up and down" to mix.
20. Place Spin Plate onto a new 0.5 ml Collection Plate.



21. Load approximately 650 μ l into each well of the Spin Plate and apply Centrifuge Tape.
22. Centrifuge at room temperature for 3 minutes at 4500 x g. Discard the flow through and place the Spin Plate back on the same 0.5 ml Collection Plate. Discard the Centrifuge Tape.
23. Repeat steps 20-21 until all the supernatant has been processed. Discard the final flow through.

What's happening: DNA is selectively bound to the silica membrane in the Spin Plate device in the high salt solution. Contaminants pass through the filter membrane, leaving only DNA bound to the membrane.

24. Place the Spin Plate back on the same 0.5 ml Collection Plate.
25. **Confirm that ethanol has been added to Solution C5-D (see step 1).** Add 500 μ l of Solution C5-D to each well of the Spin Plate. Apply Centrifuge Tape to the Spin Plate.

What's happening: Solution C5-D is an ethanol based wash solution used to further clean the DNA that is bound to the silica membrane in the Spin Plate. This wash solution removes residues of salt, humic acid, and other contaminants while allowing the DNA to stay bound to the silica membrane. This flow through is waste containing ethanol wash solution and contaminants that did not bind to the silica Spin Plate membrane.

Note: You can wash more than one time to further clean DNA if desired. In some cases where soils have very high humic acid content, it will be beneficial to repeat this wash step. There is 10% extra Solution C5-D in the bottle for this purpose. Solution C5-D is also sold separately (MO BIO Catalog# 12955-4-5D & 12955-12-5D).

26. Centrifuge at room temperature for 3 minutes at 4500 x g. Discard the flow through and place the same 0.5 ml Collection Plate beneath the Spin Plate.

What's happening: This step removes residual Solution C5-D (ethanol wash solution). It is critical to remove all traces of wash solution because it can interfere with downstream DNA applications.

27. Centrifuge again at room temperature for 3 minutes at 4500 x g. Discard the flow through.

What's happening: Once again it is important to avoid any traces of the ethanol based wash solution.

28. Carefully place the Spin Plate onto a Microplate. Remove Centrifuge Tape and discard. Allow to air dry for 10 minutes at room temperature.

What's happening: This step removes residual Solution C5-D (ethanol wash solution). It is critical to remove all traces of wash solution because it can interfere with downstream DNA applications.

29. Add 100 μ l of Solution C6 to the center of each well of the Spin Plate and apply Centrifuge Tape.

What's happening: Placing the Solution C6 (sterile elution buffer) in the center of the small white membrane will make sure the entire membrane is wetted. This will result in more efficient and complete release of the DNA from the silica Spin Plate membrane.

30. Centrifuge at room temperature for 3 minutes at 4500 x g. Remove Centrifuge Tape and discard.

What's happening: As the Solution C6 (elution buffer) passes through the silica membrane, DNA that was bound in the presence of high salt is selectively released by Solution C6 (10 mM Tris pH 8.0) which is a low salt solution.



31. Cover wells of Microplate with the Elution Sealing Mat provided. DNA is now ready for any downstream application. No further steps are required.

Prolonged storage at 4°C will result in the evaporation of eluted DNA. We recommend storing DNA frozen (-20°C or -80°C). Solution C6 does not contain EDTA. To concentrate the DNA see the Hints and Troubleshooting Guide.

Thank you for choosing the PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit.

Hints and Troubleshooting Guide

Amount of Soil to Process

This kit is designed to process 0.25 gram of soil. Recommended starting amounts for different soil types are described in the table below. For wet soils, see information under “Wet Soil Sample” below.

Soil type	Recommended maximum to process
Dry sandy soil	0.5 gm
Dry clay	0.25 gm
Wet clay	0.25 gm
Potting soil	0.1 gm
Sediment	0.25 gm
Loam	0.25 gm
Peat moss	0.1 gm
Rich farm soil	0.25 gm
Amended soil	0.150 gm
Compost samples	0.1 gm

For larger sample sizes, up to 5 grams, try using our single prep PowerMax[®] Soil DNA Isolation Kit (MO BIO Catalog# 12988-10).

Wet Soil Sample

If soil sample is high in water content, weigh the slurry and dispense into the wells. It is suggested to restrict the starting amount to 0.25 g. Increasing the amount used will increase the amount of volume in the subsequent steps.

Cells are Difficult to Lyse

When working with organisms that are proven to be difficult to lyse using mechanical or chemical methods, a 10 minute incubation at 70°C, after adding Solution C1, can be performed. Continue by proceeding with the mechanical lysis step using the 96 Well Plate Shaker.

Preparation Methods for Processing Samples (Conventional Freezing)

Add the samples to the PowerSoil[®]-htp Bead Plate and maintain at either -70°C or at -20°C until the samples are completely frozen. Immediately float the PowerSoil[®]-htp Bead Plate in a 65°C water bath. Repeat the freeze-thaw a second time and proceed with the addition of Solution C1.

Optional: After the second freeze-thaw, Solution C1 can be added along with Proteinase-K (MO BIO Catalog# 1222-2) to improve the lysis efficiency for some organisms.

Vacuum Hints

If a vacuum step seems to be taking a long time, turn off the vacuum source and lift the filter plate off the vacuum to release any back pressure. Replace the filter plate and turn the vacuum source back on. Be sure there are no air leaks around the plate. If slow vacuum continues, you can centrifuge the filter plate as an alternative. Be sure any unused wells are covered with the Sealing Tape provided.

Centrifuge with a Maximum Speed Less Than 4500 x g

Multiply the protocol time and speed to determine total x g. Divide the total by the maximum speed of your centrifuge (round up if necessary). This will be the number of minutes your centrifuge will need to run to achieve the appropriate overall force. **Example:** 10 minutes at 4500 x g = 45000. If your centrifuge has a maximum speed of 2500 x g, divide 45000 ÷ 2500 = 18 minutes of centrifugation.



Hints and Troubleshooting Guide cont.

If DNA does not PCR amplify

- Check DNA yield by gel electrophoresis and spectrophotometer reading. Template is typically added to 10 ng per reaction, although more or less may be needed depending on the reaction conditions, enzyme activity, and copy number of the target sequence.
- If DNA does not amplify after altering the amount of template in the reaction, PCR optimization (i.e. changing reaction conditions, validating primers, or testing a different polymerase) should be attempted.

Eluted DNA Sample has Color

If you are working with a soil sample that has an extremely high humic acid content, there is the unlikely possibility that your eluted solution may contain some color. If this occurs, increase the number of washes with Solution C5-D to two or three washes. If you are still encountering coloring, please contact MO BIO Laboratories, Inc. for technical assistance.

DNA Floats Out of Well When Loaded on a Gel

This usually occurs because some residual Solution C5-D remains in the final sample. Prevent this by being careful not to transfer liquid onto the bottom of the spin plate. Ethanol precipitation (described in "Concentrating the DNA") is the best way to remove residual Solution C5-D.

Concentrating the DNA

The final volume of eluted DNA will be 100 μ l. The DNA may be concentrated by adding 5 μ l of 5M NaCl and inverting 3-5 times to mix. Next, add 200 μ l of 100% cold ethanol and invert 3-5 times to mix. Incubate at -20°C for at least 10 minutes to overnight. Centrifuge at 13,000 x g for 15 minutes. Decant all liquid. Wash the DNA pellet with 70% cold ethanol. Centrifuge at 13,000 x g for 10 minutes to re-pellet the sample. Decant ethanol and dry in a speed vacuum, dessicator, or ambient air. Resuspend precipitated DNA in sterile water or sterile 10 mM Tris.

Note: This procedure must be done individually after transferring the eluted sample to a microcentrifuge tube.

Storing DNA

DNA is eluted in Solution C6 (10 mM Tris). Store the DNA at -20°C to prevent degradation. DNA can be eluted in TE without DNA loss, but the EDTA may inhibit downstream reactions such as PCR and automated sequencing. DNA may also be eluted with sterile DNA-Free PCR Grade Water (MO BIO Catalog# 17000-10). DNA that has been eluted into sterile water should be stored at -70°C. Prolonged storage in the Microplates at 4°C will result in the loss of liquid due to evaporation.

MO BIO offers TE-4 (10 mM Tris, 0.1 mM EDTA pH 8.0) which will allow for maximal protection of DNA during storage with no PCR inhibition (Catalog# 17320-1000).



Additional Applications and References

Root tips/Mycorrhizae

Microbial colonization of root tips and microbial growth penetrating the root tips, are both common phenomenon occurring in many plant species. For optimal processing of these samples, please follow the suggested sample processing method above. Transfer the powdered mass into the PowerSoil[®]-htp Bead Plate supplied with this kit and proceed with the soil protocol. This procedure ensures that organisms colonizing the root tips, both externally and internally, are released for efficient lysis.

Reference

De Souza, F.A., G.A. Kowalchuk, P. Leeflang, J.A. van Veen, and E. Smit. 2004. PCR-Denaturing Gradient Gel Electrophoresis Profiling of Inter- and Intraspecies 18s rRNA Gene Sequence heterogeneity is an Accurate and Sensitive method to Assess Species Diversity of Arbuscular Mycorrhizal Fungi of the Genus *Gigaspora*. *Appl. Environ. Microbiol.* 70 (3): 1413-1424.

Microbial Mats and Biofilms

Microbial mats and biofilms are a common occurrence in many natural environments. The mats or biofilms provide protection for the microbial community that is growing within the biofilms; direct lysis of the microbial cells within this protected community is inefficient. For processing biofilm and biomat samples, we recommend the PowerBioFilm™ DNA and RNA kits. For more information go to www.mobio.com/biofilm.

Fungal Mats

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit is suitable for isolating DNA from fungi grown in pure culture. Harvesting a freshly grown culture is suggested because the growing hyphal tips of the fungal mat are a rich source of DNA. This is accomplished by harvesting the fungal mat using cheesecloth; this will separate single cell organisms from the fungal mat. After filtering, transfer an appropriate amount of the retained moist fungal mat into the PowerSoil[®]-htp Plate. To improve the lysing efficiency of this kit, freeze/thaw cycles may be performed after adding samples to the bead plate and adding the Bead Solution. Store the plate at -70°C or -20°C until the samples are completely frozen. Immediately float the PowerSoil[®]-htp Bead Plate in a 65°C water bath to thaw. Repeat the freeze-thaw a second time and proceed with the addition of Solution C1.

Spores (bacterial and fungal) in Soils

Fungal spores are tougher than bacterial spores because of their high polysaccharide content. Follow the freeze-thaw protocol detailed above and heating the samples to 70°C for 15 minutes before bead beating may also be helpful. When dealing with spores harvested from pure cultures, please use the UltraClean[®] Microbial DNA Isolation Kit (MO BIO Catalog# 12224-50).

Fecal Samples and Cattle Manure

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit can be used without significant modifications on a wide variety of fecal materials.

References

1. Clement, B.G., and C.L. Kitts. 2000. Isolating PCR Quality DNA from Human Feces with a Soil DNA Kit. *Biotechniques* 28(4): 640-644.
2. Ibekwe, A.M., P. Watt, C. Grieve, V. Sharma, and S. Lyons. 2002. Multiplex Fluoregenic Real-Time PCR for Detection and Quantification of *Escherichia coli* O157:H7 in Dairy Wastewater Wetlands. *Appl. Environ. Microbiol.* 68(10): 4853-4862.
3. Ibekwe, A.M., C.M. Grieve, and S.R. Lyon. 2003. Characterization of Microbial Communities and Composition in Constructed Dairy Wetland Wastewater Effluent. *Appl. Environ. Microbiol.* 69(9): 5060-5069.
4. Trochimchuk, T., J. Fotheringham, E. Topp, H. Schraft, and K.T. Leung. 2003. A comparison of DNA extraction and purification methods to detect *Escherichia coli* O157:H7 in cattle manure. *J. Microbiol. Methods* 54: 165-175.



Additional Applications and References cont.

Rumen Samples

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit can be used to extract PCR-quality DNA from rumen samples as described in the reference below. Rumen has an actively metabolizing microbial community; unfortunately it is also rich in phenolic compounds and polysaccharides. These contaminants often inhibit downstream molecular analyses. Using the MO BIO Laboratories UltraClean[®] Soil DNA Isolation Kit or PowerSoil[®] DNA Isolation Kit, with the modifications in the publication below, will result in application ready DNA.

Reference

Mackie, R.I., R.I. Aminov, W. Hu, A.V. Klieve, D. Ouwerkerk, M.A. Sundest, and Y. Kamagata. 2003. Ecology of Uncultivated Oscillospira Species in the Rumen of Cattle, Sheep and Reindeer as Assessed by Microscopy and Molecular Approaches. Appl. Environ. Microbiol. 69(11): 6808-6815.

Insects

The UltraClean[®] Tissue & Cells DNA Isolation Kit, (MO BIO Catalog# 12334-50), is recommended for this type of starting material.

Coral Reef

The PowerSoil[®]-htp 96 Well Soil DNA Isolation Kit is also suitable for isolating PCR-quality DNA from microbial communities living within Coral reef samples. Briefly, 1.3 cm diameter core samples were drilled from coral reefs and the microorganisms were washed off the cores in 10X TE buffer. Both the slurry and pelleted microbial cells were directly used to extract microbial DNA using the UltraClean[®] Soil DNA Isolation Kit. The methodology employed is described in more detail in the following reference.

Reference

Rohwer, F., M. Breitbart, J. Jara, F. Azam, and N. Knowlton. 2001. Diversity of bacteria associated with the Caribbean coral *Montastraea franksi*. Coral Reefs 20: 85-91.

Note: Kit components have been used with the sample types listed above. This 96 Well version of the kit has been developed based on the success of the single prep kit.



Contact Information

Technical Support:

Phone MO BIO Laboratories, Inc. Toll Free 800-606-6246, or 760-929-9911

Email: technical@mobio.com

Fax: 760-929-0109

Mail: MO BIO Laboratories, Inc, 2746 Loker Ave West, Carlsbad, CA 92010

Ordering Information:

Direct: Phone MO BIO Laboratories, Inc. Toll Free 800-606-6246, or 760-929-9911

Email: orders@mobio.com

Fax: 760-929-0109

Mail: MO BIO Laboratories, Inc, 2746 Loker Ave West, Carlsbad, CA 92010

For the distributor nearest you, visit our website at www.mobio.com/distributors



Other Quality Products Available from MO BIO Laboratories, Inc.

For more product and detailed information go to www.mobio.com/catalog-request to request a catalog.

DNA Purification and Gel Extraction	Catalog No.	Quantity
PowerClean® DNA Clean-Up Kit	12877-50	50 preps
UltraClean® 15 DNA Purification Kit	12100-300	300 preps
UltraClean® PCR Clean-Up Kit	12500-50 12500-100 12500-250	50 preps 100 preps 250 preps
UltraClean®-htp 96 Well PCR Clean-Up Kit	12596-4 12596-12	4 x 96 preps 12 x 96 preps
UltraClean® GelSpin® DNA Extraction Kit	12400-50 12400-100 12400-250	50 preps 100 preps 250 preps
Plasmid DNA Isolation	Catalog No.	Quantity
UltraClean® 6 Minute Mini Plasmid Prep Kit	12300-50 12300-100 12300-250	50 preps 100 preps 250 preps
UltraClean® Standard Mini Plasmid Prep Kit	12301-50 12301-100 12301-250	50 preps 100 preps 250 preps
UltraClean®-htp 96 Well Plasmid Prep Kit	12396-4 12396-12	4 x 96 preps 12 x 96 preps
UltraClean® Midi Plasmid Prep Kit	12700-20 12700-50	20 preps 50 preps
UltraClean® Maxi Plasmid Prep Kit	12600-10 12600-20	10 preps 20 preps
UltraClean® Endotoxin-Free Mini Plasmid Prep Kit	12311-100 12311-250	100 preps 250 preps
UltraClean® Endotoxin-Free Midi Plasmid Prep Kit	12711-10	10 preps
UltraClean® Endotoxin-Free Maxi Plasmid Prep Kit	12611-10	10 preps
UltraClean® Endotoxin Removal Kit	12615	1 kit
UltraClean® Endotoxin-Free Ethanol Precipitation Kit	12616	1 kit
UltraClean® Endotoxin Removal Reagent	12625-25	25 ml
Endotoxin-Free Sodium Chloride	12626-15	15 ml
Endotoxin-Free Centrifuge Tubes	12617-100 12618-50 12619-25	100 each/2 ml tubes 50 each/15 ml tubes 25 each/50 ml tubes
RNA Isolation	Catalog No.	Quantity
PowerWater® RNA Isolation Kit (No Filters)	14700-50-NF	50 preps
PowerLyzer™ UltraClean® Tissue & Cells RNA Isolation Kit	15055-50	50 preps
PowerLyzer™ UltraClean® Plant RNA Isolation Kit	13355-50	50 preps
PowerBiofilm™ RNA Isolation Kit	25000-50	50 preps
LifeGuard™ Soil Stabilization Solution	12868-100 12868-1000	100 ml 1 L
On-Spin Column DNase I Kit (RNase-Free)	15100-50	50 preps
BiOstic® Stabilized Blood RNA Isolation Kit	12231-50 12231-100	50 preps 100 preps
BiOstic® Blood Total RNA Isolation Kit	12230-50	50 preps

RNA Isolation ... Continued	Catalog No.	Quantity
RNA PowerSoil® DNA Elution Accessory Kit	12867-25	25 preps
RNA PowerSoil® Total RNA Isolation Kit	12866-25	25 preps
UltraClean® Microbial RNA Isolation Kit	15800-50 15800-250	50 preps 250 preps
UltraClean® Tissue & Cells RNA Isolation Kit	15000-50 15000-250	50 preps 250 preps
UltraClean® Plant RNA Isolation Kit	13300-20 13300-50	20 preps 50 preps
Genomic DNA Isolation	Catalog No.	Quantity
PowerLyzer™ PowerSoil® DNA Isolation Kit	12855-50 12855-100	50 preps 100 preps
PowerLyzer™ UltraClean® Microbial DNA Isolation Kit	12255-50	50 preps
PowerBiofilm™ DNA Isolation Kit	24000-50	50 preps
PowerFood™ Microbial DNA Isolation Kit	21000-50 21000-100	50 preps 100 preps
BiOstic® Bacteremia DNA Isolation Kit	12240-50	50 preps
BiOstic® FFPE Tissue DNA Isolation Kit	12250-50	50 preps
BiOstic® Paraffin Removal Reagent	12251-50	2 x 25 ml
PowerMax® Soil DNA Isolation Kit	12988-10	10 preps
PowerSoil® DNA Isolation Kit	12888-50 12888-100	50 preps 100 preps
PowerSoil®-htp 96 Well Soil DNA Isolation Kit	12955-4 12955-12	4 x 96 preps 12 x 96 preps
UltraClean® Soil DNA Isolation Kit	12800-50 12800-100	50 preps 100 preps
UltraClean®-htp 96 Well Soil DNA Isolation Kit	12896-4 12896-12	4 x 96 preps 12 x 96 preps
UltraClean® Mega Soil DNA Isolation Kit	12900-10	10 preps
PowerClean® DNA Clean-Up Kit	12877-50	50 preps
UltraClean® Fecal DNA Isolation Kit	12811-50 12811-100	50 preps 100 preps
PowerMicrobial® Midi DNA Isolation Kit	12225-25	25 preps
PowerMicrobial® Maxi DNA Isolation Kit	12226-25	25 preps
UltraClean® Microbial DNA Isolation Kit	12224-50 12224-250	50 preps 250 preps
UltraClean®-htp 96 Well Microbial DNA Isolation Kit	10196-4 10196-12	4 x 96 preps 12 x 96 preps
PowerPlant® DNA Isolation Kit	13200-50 13200-100	50 preps 100 preps
UltraClean® Plant DNA Isolation Kit	13000-50 13000-250	50 preps 250 preps



Other Quality Products Available from MO BIO Laboratories, Inc.

For more product and detailed information go to www.mobio.com/catalog-request to request a catalog.

Genomic DNA Isolation ...Continued	Catalog No.	Quantity
UltraClean®-htp 96 Well Plant DNA Isolation Kit	13096-4 13096-12	4 x 96 preps 12 x 96 preps
UltraClean® Tissue & Cells DNA Isolation Kit	12334-50 12334-250	50 preps 250 preps
UltraClean®-htp 96 Well Tissue DNA Isolation Kit	12996-4 12996-12	4 x 96 preps 12 x 96 preps
UltraClean® Blood DNA Isolation Kit (Non-Spin)	12000-100	100 preps
UltraClean® Blood DNA Isolation Kit (Processes 1,000 ml of Blood)	12000-1000	1 kit
UltraClean® Blood DNA Isolation Kit Plus RNase (Processes 1,000 ml of Blood)	12002-1000	1 kit
UltraClean® BloodSpin® DNA Isolation Kit	12200-50 12200-250	50 preps 250 preps
UltraClean®-htp 96 Well BloodSpin® DNA Isolation Kit	12296-4 12296-12	4 x 96 preps 12 x 96 preps
UltraClean® Forensic DNA Isolation Kit	14000-10 14000-20	10 isolations 20 isolations
PowerWater® Sterivex DNA Isolation Kit	14600-50-NF	50 preps
PowerWater® DNA Isolation Kit (No Filters)	14900-50-NF 14900-100-NF	50 preps 100 preps
RapidWater™ DNA Isolation Kit (No Filters)	14810-50-NF 14810-100-NF	50 preps 100 preps
UltraClean® Water DNA Isolation Kit (No filters)	14800-10-NF 14800-25-NF	10 preps 25 preps
Microbiological Culture Media	Catalog No.	Quantity
TB DRY® Powder Growth Media	12105-05 12105-1 12105-5	500 g 1 kg 5 kg
LB Broth Powder Growth Media, pH 7	12106-05 12106-1 12106-5	500 g 1 kg 5 kg
LB Agar Powder Growth Media, pH 7	12107-05 12107-1 12107-5	500 g 1 kg 5 kg
LB Broth (Lennox) Powder Growth Media, pH 7	12108-05 12108-1 12108-5	500 g 1 kg 5 kg
LB Agar (Lennox) Powder Growth Media, pH 7	12109-05 12109-1 12109-5	500 g 1 kg 5 kg
Soybean-Casein Digest Medium (TSB), USP	12114-05 12114-1 12114-5	500 g 1 kg 5 kg
Soybean-Casein Digest Agar Medium (TSA), USP	12115-05 12115-1 12115-5	500 g 1 kg 5 kg
Yeast Extract	12110-05 12110-1 12110-5	500 g 1 kg 5 kg
Tryptone	12111-05 12111-1 12111-5	500 g 1 kg 5 kg
Agar, Bacteriological Grade	12112-05 12112-1 12112-5	500 g 1 kg 5 kg

Other Reagents and Lab Accessories	Catalog No.	Quantity
20 bp DNA Ladder	17020-40	40 µg
100 bp DNA Ladder	17100-40	40 µg
1 kb DNA Ladder	17200-100	100 µg
UltraClean® Agarose, Molecular Biology Grade	15003-50 15003-100 15003-500 15003-1000	50 g 100 g 500 g 1 kg
UltraClean® MS-8 Agarose	15515-50 15515-100 15515-500	50 g 100 g 500 g
UltraClean® Forensic Agarose	15505-50 15505-100 15505-500	50 g 100 g 500 g
UltraClean® Low Melt Agarose	15005-50 15005-100 15005-500	50 g 100 g 500 g
UltraClean® Low Melt Sieve Agarose	15004-50 15004-100 15004-500	50 g 100 g 500 g
Ethidium Bromide Solution	15006-1 15006-10	1 ml 10 ml
Ethidium Bromide Destaining Tea Bags	15007-25	25 bags
Dye Dots™ Dry Gel Loading Dye with Bromophenol Blue	15020-10 15020-20	10 plates 20 plates
Bromophenol Blue Gel Loading Buffer	15008-1 15008-5	1 ml 5 x 1 ml
Bromophenol Blue/Xylene Cyanol Gel Loading Buffer	15009-1 15009-5	1 ml 5 x 1 ml
TAE Buffer, 50X (Tris-acetate-EDTA)	15001-500 15001-1000	500 ml 1 liter
TBE Buffer, 10X (Tris-borate-EDTA)	15002-500 15002-1000	500 ml 1 liter
RNase-Free Gloves	1556-XS 1556-S 1556-M 1556-L	bag of 150 bag of 150 bag of 150 bag of 150
UltraClean® Lab Cleaner	12095-250 12095-500 12095-1000	250 ml squeeze bottle 500 ml spray bottle 1 liter bottle
KAPA PROBE FAST qPCR Kits	51220-100 51220-500 51220-1000	100 reactions 500 reactions 1000 reactions
KAPA SYBR® FAST Universal 2X qPCR Master Mix	51230-100 51230-500 51230-1000	100 reactions 500 reactions 1000 reactions
KAPA2G Robust HotStart ReadyMix	51240-100 51240-500	100 reactions 500 reactions
KAPA HiFi HotStart ReadyMix	51250-100 51250-500	100 reactions 500 reactions
KAPA2G FAST HotStart DNA Polymerase with dNTPs	51260-100 51260-250 51260-500	100 reactions 250 reactions 500 reactions



Other Quality Products Available from MO BIO Laboratories, Inc.

For more product and detailed information go to www.mobio.com/catalog-request to request a catalog.

Other Reagents and Lab Accessories... Continued	Catalog No.	Quantity
KAPA2G FAST HotStart ReadyMix	51270-100 51270-500	100 reactions 500 reactions
KAPA Long Range HotStart DNA Polymerase with dNTPs	51280-100 51280-250 51280-500	100 reactions 250 reactions 500 reactions
KAPA Taq Polymerase ReadyMix	51290-250	250 reactions
DNase (RNase-Free)	15600-5 15601-100	5 mg 2500 units
Proteinase K	1223-100 1222-2	100 mg 2 ml (20 mg/ml)
Ribonuclease A (25 mg/ml)	1202-1 1202-5	1 ml 5 ml
PCR Water	17000-1 17000-5 17000-10 17000-11	1 ml 5 x 1 ml 10 x 1 ml 10 ml bottle
Molecular Biology Grade Water	17012-200 17012-5200	200 ml 5 x 200 ml
DEPC Treated Water	17011-200 17011-5200	200 ml 5 x 200 ml
Endotoxin-Free Water	17013-10 17013-50 17013-100 17013-500	10 ml 50 ml 100 ml 500 ml
UltraClean® 5N NaOH	17300-30	30 ml
UltraClean® 0.1N NaOH	17301-30	30 ml
UltraClean® 0.5M EDTA, pH 8.0	17310-100	100 ml
UltraClean® TE-4 Buffer	17320-1000	1L
UltraClean® TE Buffer	17325-1000	1L
UltraClean® 1X PBS, pH 7.4	17330-500	500 ml
UltraClean® 2M Tris, pH 8.0	17370-250	250 ml
UltraClean® 0.1M Tris, pH 7.0	17371-1000	1L
Instrumentation and Accessories	Catalog No.	Quantity
PowerLyzer™ 24 Bench Top Bead-Based Homogenizer (110/220V)	13155	1 unit
PowerLyzer™ Tube Holder	13156	1 unit
PowerLyzer™ Tube Holder Stand	13157	1 unit
PowerVac™ Mini System	11992	1 unit + 20 adapters
PowerVac™ Manifold	11991	1 unit
PowerVac™ Mini Spin Filter Adapters	11992-10 11992-20	10 adapters 20 adapters
Ceramic Bead Tubes, 1.4 mm	13113-50	50 bead tubes
Ceramic Bead Tubes, 2.8 mm	13114-50	50 bead tubes
Glass Bead Tubes, 0.5 mm	13116-50	50 bead tubes
Glass Bead Tubes, 0.1 mm	13118-50	50 bead tubes
Metal Bead Tubes, 2.38 mm	13117-50	50 bead tubes
2.0 ml Tough Tubes with Cap	13119-500 13119-1000	500 1000
Carbide Bead Tubes, 0.25 mm	13121-50	50 x 0.5 ml tubes
Garnet Bead Tubes, 0.15 mm	13122-50	50 x 0.5 ml tubes

Instrumentation and Accessories... Continued	Catalog No.	Quantity
Garnet Bead Tubes, 0.70 mm	13123-50	50 x 2 ml tubes
Garnet + ¼ Ceramic 15 ml Bead Tubes, 0.70 mm	13134-50	50 tubes
Garnet + ¼ Ceramic 50 ml Bead Tubes, 0.70 mm	13144-10 13144-50 13144-100 13144-500	10 tubes 50 tubes 100 tubes 500 tubes
Glass 15 ml Bead Tubes, 0.1 mm	13135-50	50 tubes
Glass 50 ml Bead Tubes, 0.1 mm	13145-10 13145-50 13145-100 13145-500	10 tubes 50 tubes 100 tubes 500 tubes
Glass 15 ml Bead Tubes, 1.0 mm	13136-50	50 tubes
Ceramic 15 ml Bead Tubes, 1.4 mm	13137-50	50 tubes
Ceramic 50 ml Bead Tubes, 1.4 mm	13147-10 13147-50	10 tubes 50 tubes
Metal 50 ml Bead Tubes, 2.38 mm	13149-10 13149-50	10 tubes 50 tubes
Ceramic Beads, 1.4 mm	13113-325	325 gm
Ceramic Beads, 2.8 mm	13114-325	325 gm
Glass Beads, 0.5 mm	13116-400	400 gm
Metal Beads, 2.38 mm	13117-500	500 gm
Glass Beads, 0.1 mm	13118-400	400 gm
Carbide Beads, 0.25 mm	13121-500	500 gm
Garnet Beads, 0.15 mm	13122-500	500 gm
Garnet Beads, 0.70 mm	13123-05	500 gm
PowerMix 15 ml Bead Tubes	13138-50	50 tubes
PowerMix 50 ml Bead Tubes	13148-10 13148-50	10 tubes 50 tubes
2 ml Collection Tubes	1200-100-T 1200-150-T 1200-250-T	100 tubes 150 tubes 250 tubes
2 ml Screw Cap Tubes	12800-200-E	200 tubes & caps
15 ml Collection Tubes	12700-T	25 tubes
50 ml Centrifuge Tubes	12600-T	25 tubes
Spin Filters (in 1.9 ml tubes)	1200-50-SF 1200-100-SF 1200-250-SF	50 filters 100 filters 250 filters
Endotoxin-Free Centrifuge Tubes	12617-100 12618-50 12619-25	100 each/2 ml tubes 50 each/15 ml tubes 25 each/50 ml tubes
15 ml Midi Spin Filters	12700-SF	25 spin filters



Other Quality Products Available from MO BIO Laboratories, Inc.

For more product and detailed information go to www.mobio.com/catalog-request to request a catalog.

Instrumentation and Accessories... <i>Continued</i>	Catalog No.	Quantity
Vortex-Genie® 2 Vortex (120V)	13111-V	1 unit
Vortex-Genie® 2 Vortex (220V)	13111-V-220	1 unit
Vortex Adapter, holds 12 (1.5-2.0 ml) tubes	13000-V1	1 unit
Vortex Adapter, holds 6 (5 ml) tubes	13000-V1-5	1 unit
Vortex Adapter, holds 4 (15 ml) tubes	13000-V1-15	1 unit
Vortex Adapter, holds 2 (50 ml) tubes	13000-V1-50	1 unit
Vortex Adapter, holds 24 (1.5-2.0 ml) tubes	13000-V1-24	1 unit
Anti-Static Funnels, Micro	23301-96	Pack of 96
Anti-Static Funnels, Small	23302-50	Pack of 50
Anti-Static Funnels, Medium	23303-50	Pack of 50
Anti-Static Funnels, Large	23304-20	Pack of 20
Mini Horizontal Gel System	16001	1 each
Mini Horizontal Gel Caster, 3 place	16003	1 each
Mini Horizontal Gel Tray	16004	1 each
Polycarbonate Single-sided Comb	16005 16006 16007 16008	1 mm x 3 well 1 mm x 8 well 1 mm x 10 well 1 mm x 12 well
Polycarbonate Dual-sided Comb	16013 16014 16015 16016	1 mm x 8 well/16 well 1 mm x 10 well/14 well 2 mm x 8 well/16 well 2 mm x 10 well/14 well
Teflon Single-sided Comb	16009 16010 16011 16012	1 mm x 3 well 1 mm x 8 well 1 mm x 10 well 1 mm x 12 well
Teflon Dual-sided Comb	16017 16018 16019 16020	1 mm x 8 well/16 well 1 mm x 10 well/14 well 2 mm x 8 well/16 well 2 mm x 10 well/14 well
Power Supply w/Timer, (120V)	16023	1 unit
96 Well Plate Shaker (120V)	11996	1 unit
96 Well Plate Shaker (220V)	11996-220	1 unit
Plate Adapter Set	11999	1 set
Vacuum Pump (120V)	11998	1 unit
Vacuum Pump (220V)	11998-220	1 unit